

## Improvements in Characterization of FIB Prepared Surfaces of Aluminum Using Xe<sup>+</sup> Plasma FIB

Brandon Van Leer<sup>1</sup> and Rick Passey<sup>1</sup>

<sup>1</sup> FEI Company, 5350 NW Dawson Creek Drive, Hillsboro, OR, USA


Critical to understanding the microstructure of aluminum and aluminum alloys is the careful surface preparation either by mechanical polishing or by ion beam techniques. Cross-section and S/TEM sample preparation of aluminium and aluminium alloys to characterize grain boundaries by focused ion beam (FIB) continues to be a major interest in metallurgical analysis because of FIB's ability to prepare site specific specimens while eliminating damage from mechanical polishing or electro-polishing [1]. However, gallium is particularly mobile in aluminium and known to diffuse along grain boundaries or interfaces, which can lead to embrittlement [2,3]. Often subsequent polishing in a broad beam polisher is a necessary step to improve the quality of the prepared surface for EBSD analysis. Also, characterizing large grain aluminum samples by Ga<sup>+</sup> FIB requires longer processing time when investigating multiple grains. Recent instrumentation using plasma FIB (PFIB) technology and Xe<sup>+</sup> ions offers an alternative ion species for FIB milling and increased milling rates because of its ability to deliver 30 – 40 times more current compared to Ga<sup>+</sup> FIBs. While the measured sputter rate of aluminum using Ga<sup>+</sup> and Xe<sup>+</sup> differs by about 25% (0.31 μm<sup>3</sup>/nC [Ga] and 0.41 μm<sup>3</sup>/nC [Xe]), the ability to prepare gallium free thin sections for S/TEM analysis or large area cross-sections offers a solution to FIB milling artifacts seen with a Ga<sup>+</sup> FIB as well as improving statistical analysis of grain populations.

Conventional cross-sections by FIB and S/TEM specimens of commercial grade 6061 T6 aluminum were prepared using the Helios G4 UX DualBeam™ using 30 kV Ga<sup>+</sup> ions and a Helios PFIB DualBeam using 30 kV Xe<sup>+</sup> ions. Sample quality of the cross-sections were evaluated by comparing EBSD index rates and band contrast. Sample quality of the S/TEM specimens were evaluated by measuring FIB sidewall damage and observation of Ga contamination at grain boundaries in a Themis Z™ TEM operating at 300 keV.


A comparison of FIB sidewall damage between specimens prepared in Ga<sup>+</sup> FIB and Xe<sup>+</sup> PFIB revealed a decrease in sidewall by as much as 30% depending on the accelerating voltage applied. Figure 1 shows a SEM image the finished TEM lamellae of 6061 T6 aluminum as prepared in the Helios PFIB for FIB sidewall damage. Figure 2 shows a SEM image of a finished thin section of 6061 T6 aluminum across a grain boundary. EBSD results revealed that cross-section surfaces prepared with 30 kV Xe<sup>+</sup> PFIB had better band contrast results and improved indexing rates than cross-sections prepared with 30 kV Ga<sup>+</sup> FIB.

### References:

- [1] L.A. Giannuzzi *et al.*, Mater. Res. Soc. Symp. Proc. **480** (1997) p 19.
- [2] K.A. Unocic *et al.*, Journal of Microscopy **240** (2010) p 227.
- [3] R.C. Hugo and R.G. Hoagland, Scr. Mater. **41** (1999) p 1341.

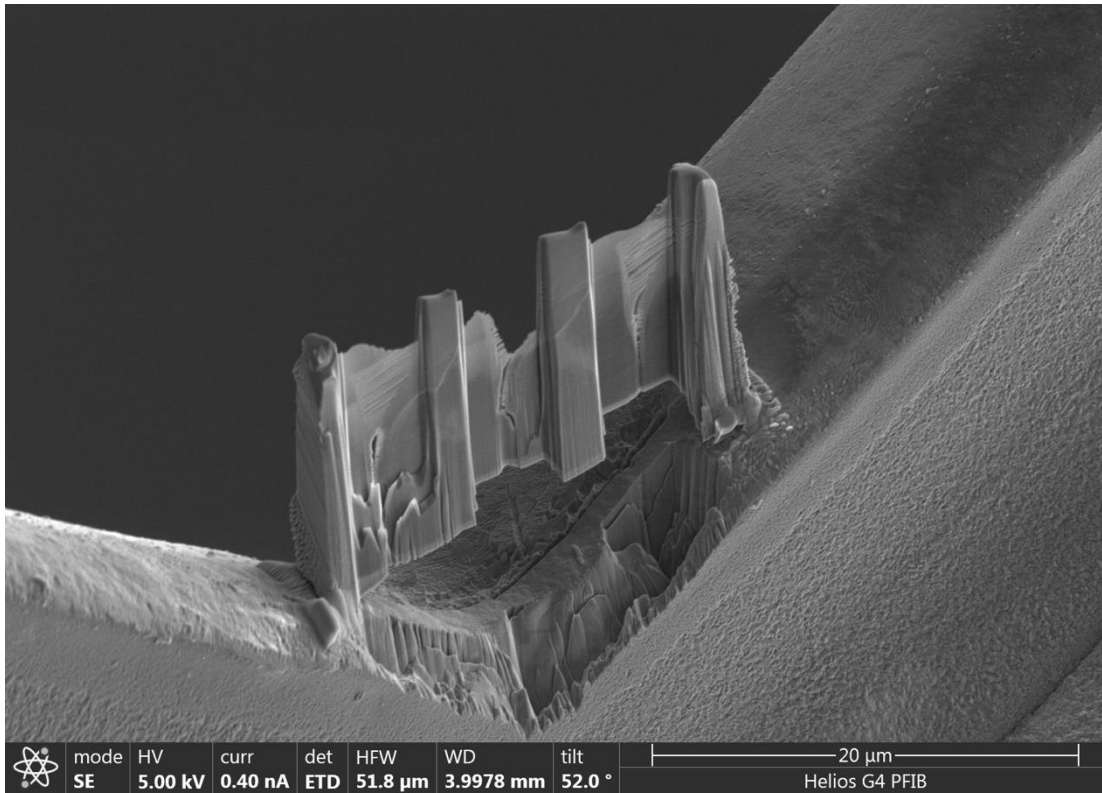

**Figure 1.** SEM image thin sectioned 6061 T6 aluminum characterizing the effect of accelerating voltage on FIB sidewall damage.

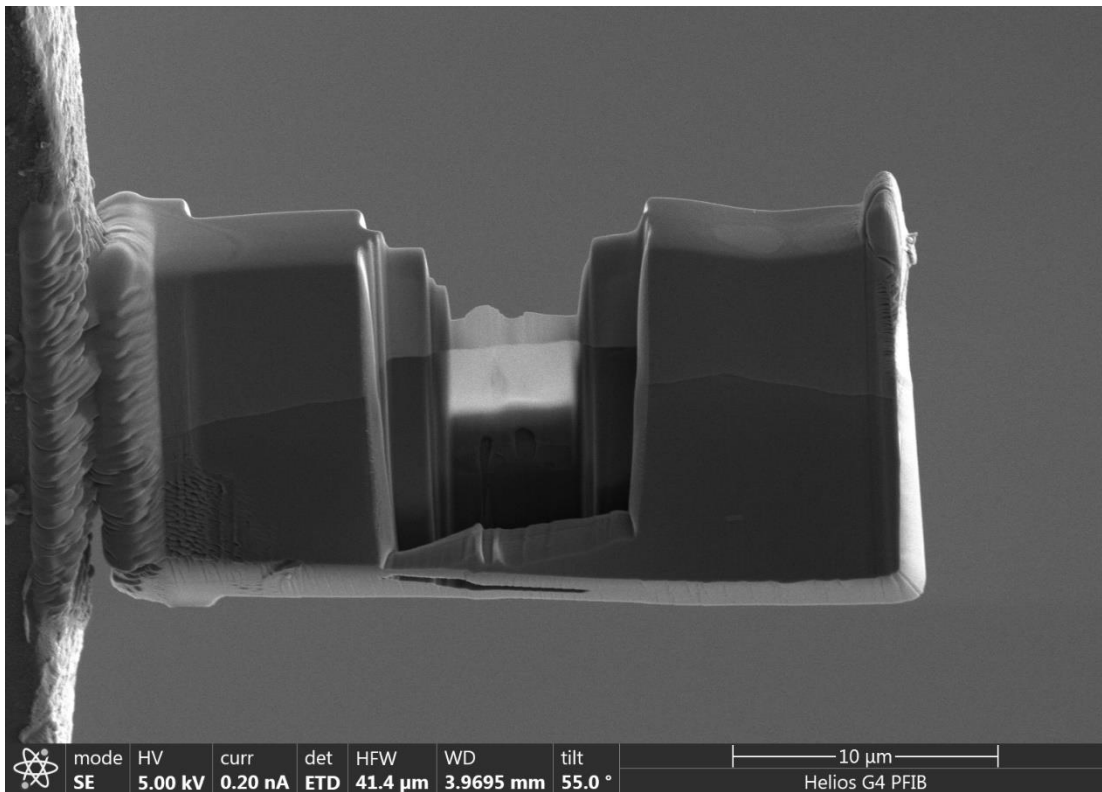

**Figure 2.** SEM image of thin sectioned 6061 T6 aluminum across a grain boundary.